- (DE) Bedienungsanleitung
- (GB) Operating Instructions
- (FR) Mode d'emploi
- (NL) Handleiding
- (IT) Istruzioni per l'uso
- (ES) Instrucciones de uso
- (PT) Manual de utilização

DE	Bedienungsanleitung	4
GB	Operating Instructions	18
FR	Mode d'emploi	30
NL	Handleiding	44
IT	Istruzioni per l'uso	58
ES	Instrucciones de uso	72
PT	Manual de utilização	86

Fig. 1

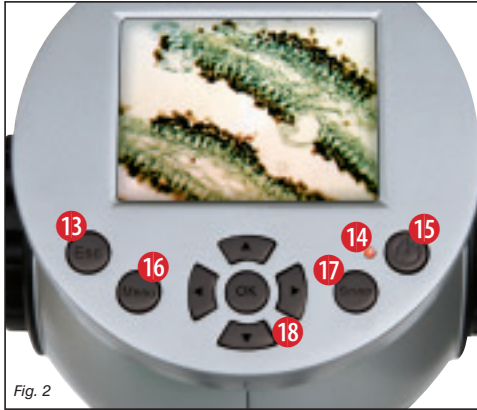
Fig. 2


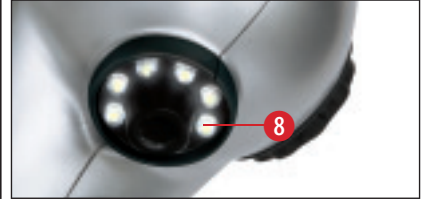
Fig. 3

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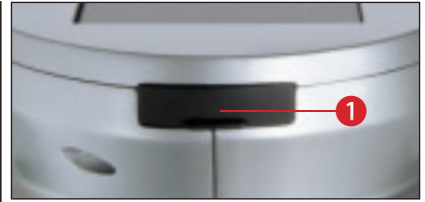RISK to your child.

 Aids with sharp edges and tips are sometimes used with this device. Please store the device and all of its accessories and aids out of the reach of children. There is a risk of **INJURY**.

This device contains electronic components that are powered by either a mains connection or batteries. Never leave a child unsupervised with this device. The device should only be used as per these instructions otherwise there is a serious **RISK** of **ELECTRICAL SHOCK**.

Batteries should be kept out of children's reach. When inserting batteries please ensure the polarity is correct. Leaking or damaged batteries can cause injury if they come into contact with the skin. If you need to handle such batteries please wear suitable safety gloves.

Children should only use this device under supervision. Keep packaging materials (plastic bags, rubber bands, etc.) away from children. There is a risk of **SUFFOCATION**.

The chemicals and liquids provided should be kept out of reach of children. Do not drink the chemicals! Hands should be washed thoroughly under running water after use.

In case of accidental contact with the eyes or mouth rinse with water. Seek medical treatment for ailments arising from contact with the chemical substances and take the chemicals with you to the doctor.

FIRE-/ DANGER OF EXPLOSION!

 Do not expose the device to high temperatures. Use only the mains adapter supplied or those battery types recommended. Never short circuit the device or batteries or throw into a fire. Exposure to high temperatures or misuse of the device can lead to short circuits, fire or even explosion!

RISK of material damage.

 Never take the device apart. Please consult your dealer if there are any defects. The dealer will contact our service centre and send the device in for repair if needed.

Do not subject the device to temperatures exceeding 60 C.

TIPS on cleaning

Remove the device from its energy source before cleaning (remove plug from socket / remove batteries)

Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Clean the lens (objective and eyepiece) only with the cloth supplied or some other soft lint-free cloth (e.g. micro-fibre). Do not use excessive pressure - this may scratch the lens.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

DISPOSAL

Dispose of the packaging material/s as legally required. Consult the local authority on the matter if necessary.

Do not dispose of electrical equipment in your ordinary refuse. The European guideline 2002/96/EU on Electronic and Electrical Equipment Waste and relevant laws applying to it require such used equipment to be separately collected and recycled in an environment-friendly manner. Empty batteries and accumulators must be disposed of separately. Information on disposing of all such equipment made after 01 June 2006 can be obtained from your local authority.

EEC conformity explanation

Meade Instruments GmbH & Co. KG has formed a “Conformity Declaration” in line with the applicable guidelines. This declaration is available for review.

These are the parts of your microscope

1. SD Card Receptacle
2. LCD Monitor
3. Objective Nosepiece
3. LCD Module
4. Mini USB Cable Receptacle
5. Objective Lens
6. Focus Knob
7. Mechanical Stage
8. surface illumination
9. substage illumination
10. Setting the brightness (-)
11. On/off key for the lighting type
12. Setting the brightness (+)
13. Esc Key
14. Monitor Control Light
15. On/off Switch (Main Switch)
16. Menu Key
17. Snap Key
18. Direction/OK Key
19. Power Adapter
20. Unit Outlet
21. Microscope Accessories
22. Software CD
23. USB Cable

1. Setting up your Microscope

Carefully remove the microscope and the other components from the package and place them on a table or other flat surface. Remove the dust cover from the microscope.

A power connection (220-230 V) is required for observation with the electric LED light and the LED eyepiece (display module).

The microscope can also connect to a PC (USB connection).

In order to operate the device, connect the supplied power plug (Fig. 3) with the power connection (Fig. 3) on the back of the microscope base (Fig. 3) to a power outlet (220-230 V). Subsequently, turn on the desired lighting using the light switch (11) and select the required brightness with the „+“ (12) and „-“ (10) keys.

2. Connection between Microscope and PC

2.1 Transferring Data to a PC

Use the USB cable to connect your PC to the microscope.

When the connection has been established successfully, the PC recognizes the new hardware on its own. The control light on the microscope turns on, but the digital camera

remains off. On your PC, open the corresponding folder for the removable device. The files for the saved pictures can be found here; copy them to your PC. You can process them further from there.

Attention:

Do not remove the USB cable during the data transfer. The data will be lost.

Before breaking the USB connection or turning off the LCD module, you need to remove the device using the hardware assistant on your PC („deactivate“).

If the drive is not deactivated, the PC could crash or you could lose your data!

2.2 Installing the Software

Attention:

Do not connect the DigiMikro DM700 to your PC (computer) yet. In order to work with the DigiMikro DM700 for the long term, we recommend that you always connect it to the same USB port. In some cases, your computer may not detect the device. If this occurs, you can typically fix the issue by re-installing the drivers (from the CD). If this does not work, please see the following chapter on troubleshooting.

Insert the supplied CD-ROM into your CD/DVD drive. The CD menu opens automati-

cally. If the menu does not open, select the CD/DVD drive in Windows Explorer. There, start the file „Menu.exe“ with a double click.

2.3 Installing the Drivers

Please make sure that the DigiMikro DM700 is not connected to your PC.

Take the follow steps:

1. Save the current data and close any open programs.
2. Click on „Install Driver“.
3. A selection screen appears. Select the language and click „Continue“.
4. A „Welcome“ screen appears. Click „Continue“.
5. The „Setup Status“ window now appears. The program files are copied to your PC. This process can take a few minutes.
6. After the window disappears, it can take a few moments before the „InstallShield Wizard Complete“ window appears. Click on „Finish“. Your PC will restart.
7. Once the computer restarts, the driver is installed.
8. Connect the supplied USB cable to a free USB port.
9. Press and hold the left arrow key (◀) on the LCD module. Now, connect the USB cable to the DigiMikro DM700. The webcam function will now start.

Driver Assignment:

The DigiMikro DM700 is now assigned to the driver. In Windows XP, a notification appears in the lower right: „The hardware is now installed and can be used.“

The DigiMikro DM700 is now operational.

2.4 Displaying and Saving Images from the DigiMikro DM700 on Your PC:

After successfully installing the DigiMikro DM700, you can use the webcam function by repeating Points 8. and 9.

2.5 Installing the Image Editing Software

Note: If you have already installed an image editing program, you do not need to go through with Point 2.4. In this way, only the driver for the DigiMikro DM700 is installed. If you are not sure, please continue with the following steps.

1. Insert the supplied CD-ROM into your CD/DVD drive. The CD menu opens automatically. If the menu does not open, select the CD/DVD drive in Windows Explorer. There, start the file „Menu.exe“ with a double click.
2. Click on „Install Software“.
3. A selection menu appears. Here, you can select the desired language. Confirm with „OK“.
4. Click „Continue“ in the „Welcome“ screen.

5. The License Agreement appears. Confirm with „Yes“.

6. The next two windows will ask you for the „Destination Directory“ and the „Program Folder.“ Confirm this with „Continue“.

7. The „Setup Status“ window appears. The program files are copied to your PC. This process can take a few minutes.

8. The „InstallShield Wizard Complete“ window appears. Click on „Finish“. The image editing software is now installed.

2.6 Troubleshooting: the device isn't recognised by the computer (USB)

1. This occurs in rare cases with some OEM Windows Vista or XP versions. The device is not defective. Updating the notebook or PC USB driver usually suffices.

2. If not, proceed as below. (The maker cannot accept any liability whatsoever in these cases¹).

Delete ghost devices. These are devices not actually currently connected to your computer. The reason for this is that Windows generates a system settings entry for each USB device (e.g USB Stick) on each port. It then seeks that/those device/s each time it's started. This slows down USB device recognition and Windows booting and causes non-recognition of USB ports.

Cleaning up the system environment and

deleting devices that are no longer in use helps here. To do so right click WORKPLACE (COMPUTER in Vista) and open the PROPERTIES menu. Then click the ADVANCED (ADVANCED SYSTEM SETTINGS in Vista) and then ENVIRONMENT VARIABLES. Choose the NEW option in ENVIRONMENT VARIABLES. Enter the following under VARIABLE NAME.

devmgr_show_nonpresent_devices

Under VARIABLE VALUE enter „1“. Confirm your input with OK and then reboot the computer. Change to the device manager. In the VIEW menu activate the SHOW DEVICES OPTION. The ghost devices previously not shown will now be displayed in light grey. Check the various categories including USB, memory size, etc. Only delete entries for devices you no longer use.¹

¹Source: Microsoft Knowledge Base
<http://support.microsoft.com/kb/315539/de>

3. Using the Microscope

3.1 Basic Prerequisites

Before you observe objects or specimens, you must turn on the LCD monitor and the illuminator. You also have to know how to operate the mechanical stage. Then you will be ready to start observing.

3.2 LCD Module

This digital microscope is different than a conventional microscope. The LCD monitor replaces the eyepieces that are used on typical microscopes for observing. As a result, it is possible to share your observations with others by using the monitor. Before you start observing, you must turn on the LCD monitor by pressing the on/off switch (see Fig. 3). Press and hold the button until the LCD monitor activates. This is the sign that the LCD is activated. Basically, this is all you have to do to get the LCD monitor ready for observing. The different keys of the LCD module are used mainly for taking pictures (photographs and videos) and they will be explained later in these instructions.

3.3 Illumination

With this device, a transmitted-light and reflected-light microscope, you are able to observe transparent as well as non-transparent objects. The image information of the given object is „transported“ by the light. The correct lighting determines whether or not you can see the object!

The Digital Microscope offers you three illumination options in total – their use depends on the object you would like to observe.

a) surface illumination (light from above),

b) substage illumination (light from below) or
c) surface and substage illumination (light from above and below at the same time).

If you press the On/Off button (11) once, the surface illumination (8) will be turned on. If you press the button a second time, the illumination switches to substage (9). If you then press the button a third time, both surface and substage illumination are engaged. Another press of the button turns the lighting completely off again.

Choosing the right lighting depends on what you would like to look at.

If you observe non-transparent (opaque objects (e.g. small animals, parts of plants, stones, coins, etc.) with this microscope, the light falls onto the object from above (reflected-light microscopy). The light then bounces back and enters the objective, which magnifies the image. The light then passes into the electronic eyepiece, which magnifies the image further.

For transparent objects (e.g. protozoa), on the other hand, the light shines from below through the opening in the stage and then through the object (transmitted-light microscopy). The light again passes through the objective and into the electronic eyepiece.

Using both lights at the same time only makes sense for semi-transparent objects. This operating mode is not recommended for trans-

parent objects on slides, since it can cause reflections.

Specimens of varying size, thickness or color require different illumination levels. There are two ways to modify the light levels during observation: the „+“ and „-“ keys on the device (press and hold the keys); or the EV function on the LCD monitor. The EV function (EV = exposure value brightness) increases or reduces the brightness of the image on the LCD screen. This is controlled by the right or left arrow key (▶, ◀) (Fig. 2).

If the specimen is not transparent or is dark in color, you typically need more light in order to display certain characteristics. To do so, it is best to increase the light intensity by pressing „+“ on the device.

You can find optimal lighting by experimenting with various settings, since every object requires different lighting.

3.4 Observing an Object

Using the focus wheel (Fig. 1, No. 6), move the stage all the way down. Start with a simple observation. For this, place a prepared slide on the stage, directly under the objective. Move the stage clips to the side, place the prepared slide on the stage and carefully place the stage clips onto the slide. The ob-

ject must be situated directly under the lighting.

Focus the image using the focus wheel (Fig. 1, No. 6).

You can increase the magnification with the digital zoom. To do so, use the up and down arrow keys (▲, ▼) on the LCD module. The current zoom setting is displayed on the screen.

4. Digital Pictures

Basic Operations

Activation

If the unit is off, push the on/off switch (see Fig. 2) and hold it for 1 or 2 seconds.

Deactivation

If the unit is on, push the on/off switch (see Fig. 2) and hold it for 1 or 2 seconds.

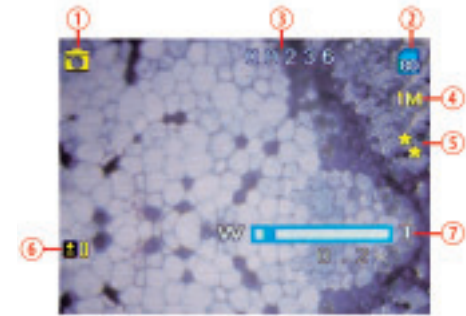
Adjustment

To change the current unit settings, call up the menu by pressing the menu key. Push the direction keys to select the corresponding settings. After choosing a setting, press the OK key to confirm it.

Display on the LCD Monitor


By pressing the Esc key once or multiple times, you access different operating modes. These are described in the following.

Digital Photo Mode


1. Digital photo mode symbol
2. Memory Symbol. If you have inserted an SD memory card, a corresponding symbol appears.
3. Number of pictures that can be stored in the memory
4. Picture resolution (number of pixels) 3M, 2M, 1M, VGA
5. Quality, ***, **, *
6. EV, exposure value brightness
7. Magnification

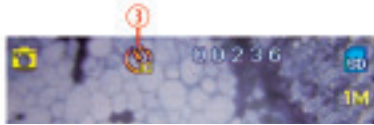Self-timer Mode


1. Self-timer mode symbol

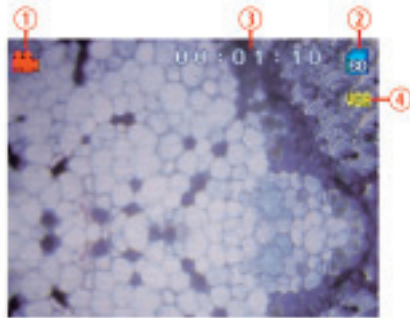Digital Video Mode

1. Digital video mode symbol
2. Memory Symbol. If you have inserted an SD memory card, a corresponding symbol appears.
3. Amount of video that can be stored in the memory
4. Picture resolution (number of pixels) VGA, QVGA

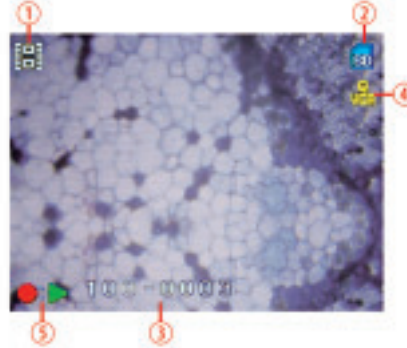Picture Index Mode

1. Symbol for photo or video
2. Memory Symbol. If you have inserted an SD memory card, a corresponding symbol appears.
3. Name of the currently selected file
4. Image resolution
5. Symbol for start, stop and pause for video mode

5. Settings

Resolution Setting

- Press the menu key in either digital photo, digital video or self-timer mode in order to open the settings menu.
- Select the required resolution (3M = 2048×1536, 2M = 1600×1200, 1M =

1280×1024, VGA = 640×480, QVGA = 320×240 for videos) with the arrow keys (▲, ▼).

- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

The resolution depends on the type of picture that you want. The higher the resolution, the better the picture quality. However, the required storage space is also higher, so fewer pictures can be saved. It is best to use the available flash memory with a carefully adjusted picture resolution.

As listed below, you have five options for the picture resolution (selection options). If you would like better image quality, select a higher resolution. The possible number of pictures depends on the circumstances (See table.)

Light Settings

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the light settings menu.
- Press the arrow keys (◀, ▶) to access the setup menu
- Using the arrow keys (▲, ▼), select the „Lighting“ option. Press the OK key to confirm.
- Using the arrow keys (▲, ▼), select the

Resolution	Storage Space				
	64M	128M	256M	512M	1G
2048×1536	120	240	480	960	1920
1600×1200	170	340	680	1360	2720
1280×1024	230	460	920	1840	3680
640×480	430	860	1720	3440	6880

Unit: one photo

desired EV value (-3, ... 0, ... +3). EV = -3 produces a very dark image; EV = +3 produces a very bright one.

- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

Focus Settings

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Sharpness“ option. Press the OK key to confirm.
- Use the arrow keys (▲, ▼) to select the required sharpness (normal, sharp, soft).
- Press „OK“ to confirm and „Snap“ to leave the menu.

- Press „Snap“ to leave the menu without changes.

Setting the Image Quality

- In the digital photo, digital video or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Quality“ option. Press the OK key to confirm.
- Use the arrow keys (▲, ▼) to select the required quality (super, fine, normal).
- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

The higher the image quality, the more drive space is required, which means fewer images

can be saved. It is best to select a reasonable image quality based on the available storage space.

Setting the Image Labeling

The device has an image-labeling function (time of exposure) for working with the images. Please set the date and time before using this function.

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Date“ option. Press the OK key to confirm.
- Using the arrow keys (▲, ▼), select either „Off“ (no image labeling) or „On“ (image labeling with date and time). Press OK to confirm.
- Using the arrow keys (▲, ▼), select the „Set Date“ option. Press the OK key to confirm. You are now in the entry menu.
- Use the left and right arrow keys (◀, ▶) to select the corresponding entry (date, time, date display) and select the setting with the up and down keys (▲, ▼).
- After the entry, press OK to confirm.

Automatic Deactivation

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Auto Off“ option. Press the OK key to confirm.
- Select the desired setting (1 minute, 2 minutes, 3 minutes, ?) with the arrow keys (▲, ▼).
- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

Language Settings

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Language“ option. Press the OK key to confirm.
- Use the arrow keys (▲, ▼) to choose one of the 8 languages.
- Press „OK“ to confirm and „Snap“ to leave the menu.

- Press „Snap“ to leave the menu without changes.

Frame Rate

In order that no image distortions appear in a video, you can select a frame rate of either 50Hz or 60Hz.

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Frequency“ option. Press the OK key to confirm.
- Use the arrow keys (▲, ▼) to select the frequency (50Hz, 60Hz).
- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

Format Function

Attention: Formatting deletes all files!

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the

„Format“ option. Press the OK key to confirm.

- Use the arrow keys (▲, ▼) to select either „Cancel“ (settings are not changed) or „OK“ (settings are reset).
- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

Default Settings

The default settings function resets all settings to their original setup.

- In the digital photo or self-timer modes, select the menu key (Menu) in order to open the settings menu.
- Press the arrow keys (◀, ▶) in order to access the setup menu
- Using the arrow keys (▲, ▼), select the „Default Settings“ option. Press the OK key to confirm.
- Use the arrow keys (▲, ▼) to select either „Cancel“ (settings are not changed) or „OK“ (settings are reset).
- Press „OK“ to confirm and „Snap“ to leave the menu.
- Press „Snap“ to leave the menu without changes.

Defaults:

Resolution	1M
Lighting	0
Sharpness	Normal
Image Quality	Fine
Date Display	YYYY/MM/DD
Auto Off	2 minutes
Language	English
Frame rate	60Hz

Functions

Digital Photo Function

Before taking digital photographs, please adjust the corresponding settings.

Attention: The number of pictures that can be taken is proportional to the free memory storage space.

- When you turn on the unit, the digital photo mode is the standard mode.
- Simply press the snap key to take a picture.

Self-timer Function

Before using the self-timer function, you need to select the corresponding settings in the self-timer menu.

- Pressing the Esc key takes you to the self-timer menu. (In digital photo mode, press-

ing the Esc key once takes you to the self-timer menu.)

- Press Snap to start taking a picture. The self-timer will run for 10 seconds before taking a picture.

Digital Video Function

Before using the digital video function, select the corresponding settings.

Attention: The amount of video that can be recorded depends on the free drive space.

- Pressing the Esc key takes you to the digital video menu. (In digital photo mode, pressing the Esc key twice takes you to the digital video menu.)
- Press Snap to start recording.
- Press Snap again to stop recording.

Digital Zoom and EV Value Settings

Sometimes, the picture of an object is too small or the illumination is inadequate in order to get a satisfactory image on the LCD monitor. Here, the EV value and digital zoom (electronic magnification) settings are very helpful.

- Press the left and right direction keys (◀, ▶) to change the EV value. The EV value range goes from -3.0 to +3.0.
- Press the up and down keys (▲, ▼) to set the electronic magnification factor. This

value range goes from 0.1 to 4.00.

6. Observing Digital Photos or Videos

- You access the picture index mode by pressing the Esc key. (You press the Esc key three times to access the picture index mode when you are in the digital photo mode).
- All the saved digital photographs and digital videos (video symbol: video camera) are displayed with the aide of the monitor.
- You can browse through the list with the direction keys (◀, ▶).
- Use the direction keys (◀, ▶) to select a digital photo or video.
- If you would like to play a digital video, select the video and press the Snap key and then the up key (▲). When you press the up key once again, the playback is interrupted.
- Pres the Esc key to go back to the digital photo mode from the picture index mode.

Functions

File Protection

The unit has a protection function to prevent the unwanted deletion of saved files.

- You access the picture index mode by pressing the Esc key. (You press the Esc key three times to access the picture index

mode when you are in the digital photo mode).

- Press the menu key to open the storage menu.
- Using the direction keys (▲ , ▼), select the option „Protection.“ Press the OK key to confirm.
- Use the direction keys (▲ , ▼) to select “Single,” “All” and „Page.“ Press OK to confirm.
- Using the direction keys (▲ , ▼), select the option „Set.“ Press the OK key to confirm.

The file protection is displayed through the lock symbol. The protection can be canceled with “Cancel” and OK for individual files (in Protection / Single) or for all files (in Protection / All).

Deleting Files

- You access the picture index mode by pressing the Esc key. (You press the Esc key three times to access the picture index mode when you are in the digital photo mode).
- Press the menu key to open the storage menu.
- Using the direction keys (▲ , ▼), select the option „Delete.“ Press the OK key to confirm.
- Use the direction keys (▲ , ▼) to select

“Single,” and “All.” Press OK to confirm.

Single: Only the file with the yellow border is deleted.

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Cropping a Picture

With the „Crop Image“ function, you can select a portion of a digital photo and save it as a new image.

- Select a digital photo with the arrow keys (◀ , ▶).
- Press the menu key to open the memory menu.
- Using the arrow keys (▲ , ▼), select the „Crop Image“ option.
- Use the arrow keys (▲ , ▼) to select from „Crop“ or „Cancel“.

- If you have selected „Crop“, a new menu appears.
- Use the arrow keys (▲ , ▼) to choose the appropriate section (green box).
- Press the Snap key to confirm your selection.
- The selection is now shown on the LCD display as a new image.
- Press the Snap key to save the image or the Menu key to exit this function without saving.

Using an SD Card

How to use the SD card

- If you have not inserted a SD card, the files will be saved in the flash memory.
- When inserting a SD card make sure the contacts are facing up. Do not push the card in forcefully. This can damage the SD card.
- The LCD monitor displays a SD card symbol. The storage space available on the SD card is displayed on the bottom right of the monitor.
- Before you remove the card, the digital microscope must be shut off.

7. Starting and Using the Photomizer SE Software

1. Start the Photomizer SE software on your computer.
2. For the best image result with the Photomizer SE software, we recommend setting the resolution at 1278x960. To do this, click on „adjust resolution“. If the resolution is not available, you can add it using the (+) field next to the resolution selection.
3. In the program window, click on „Open camera“.
4. If you have more than one device connected to your PC, you can choose the desired device in the next selection. If you only have one connected device, you do not need this step.
5. Now, you should be able to see the image from the camera on your monitor. On the microscope, bring the image into focus by turning the focus adjustment knobs (6).
6. Click on „Capture“ in order to capture an image you would like to save. This will be shown on the right in the menu bar.
7. Select this image with a single mouse click, then click on „Transfer image“.
8. You will now leave the image-capture stage, and go back to the program window.
9. Now click at the top of the screen on “File” and then on “save file” in order to secure the image.

Tip:

If you have questions or problems with the “Photomizer SE” software, you can always go to the program’s menu bar, and click on „?“ and then on „Open help“. For questions or problems, you can look on the website www.photomizer.net. Please ask your parents before using the Internet.

8. Accessories

Cover slips

You use the cover slips to cover specimens that you have placed on a slide in order to observe with a microscope.

Slides

You place the specimen on a slide. After you have covered it with some water or gum media, you can place a cover slip on top.

Tweezers

These are a type of pincers that you can use to pick up smaller objects.

Dissecting needle

A dissecting needle can be used for many things. You can use it to scratch the surface of a specimen, to fix objects in place or to stir liquids.

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