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Art. No. 9810301

- DE Bedienungsanleitung
- EN Operating instructions

## SERVICE AND WARRANTY:

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EN

## Operating Instructions

**WARNING!** Not for children under three years. Choking hazard – small parts. Contains functional sharp edges and points!

**ATTENTION:** Only suitable for children of at least 8 years of age. Instructions for parents or other responsible persons are enclosed and must be followed. Keep packaging as it contains important information.

### GENERAL WARNINGS

- **Choking hazard** – This product contains small parts that could be swallowed by children. This poses a choking hazard.
- **Risk of electric shock** – This device contains electronic components that operate via a power source (power supply and/or batteries). Only use the device as described in the manual, otherwise you run the risk of an electric shock.
- **Risk of fire/explosion** – Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- **Risk of chemical burn** – Make sure you insert the batteries correctly. Empty or damaged batteries could cause burns if they come into contact with the skin. If necessary, wear adequate gloves for protection.
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands or with different capacities. Remove the batteries from the unit if it has not been used for a long time.
- Never recharge normal, non-rechargeable batteries. This could lead to explosion during the charging process.
- Tools with sharp edges are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- Keep instructions and packaging as they contain important information.

## DISPOSAL

Dispose of the packaging materials properly, according to their type (paper, cardboard, etc). Contact your local waste disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per the Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

Empty old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 01.06.2006 from your local waste disposal service or environmental authority.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please pay attention to dispose of your used batteries as required by law - at a local collection point or in the retail market (a disposal in domestic waste violates the Battery Directive).

Batteries that contain toxins are marked with a sign and a chemical symbol. „Cd“ = cadmium, „Hg“ = mercury, „Pb“ = lead.

Cd<sup>1</sup>

Hg<sup>2</sup>

Pb<sup>3</sup>

<sup>1</sup> battery contains cadmium

<sup>2</sup> battery contains mercury

<sup>3</sup> battery contains lead

## EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address: [www.bresser.de/download/9810301/CE/9810301\\_CE.pdf](http://www.bresser.de/download/9810301/CE/9810301_CE.pdf)

---

**Your microscope is made up of the following parts (Fig. 1-4):**

- 1 WF 10x Eyepiece
- 2 45° inclined monocular head
- 3 Objective
- 4 Illumination (LED)
- 5 Microscope stage (with object plate)
- 6 Stage clips
- 7 Focus knob
- 8 On/off switch
- 9 Battery compartment
- 10 Mounting screws
- 11 Diffusion cap

**Location**

Before you start, choose an ideal location for using your microscope. It is important that you have enough light available (window, desk lamp).

**Simple Observation**

Turn the focus knob (7) to the upper stop. Now, place an object on the stage (5), directly under the objective (1). When you take a look through the eyepiece, you can see the magnified object. At this point, you still might see a slightly fuzzy picture. Adjust the image sharpness by slowly turning the focus knob (7).

**Rotatable object plate**

Test with which plate side (black or white) your object can be viewed best.

**Illumination**

For observation with the electronic light source (4) you need to insert 2 AA batteries 1.5 V, in the battery compartment (9) on the base of the microscope. The battery compartment is opened using a Phillips screwdriver. Insert the batteries with the correct polarity (+/- indication). Put the battery cover first into the small opening so that the lid fits perfectly. Now you can tighten the screw. You can not turn on the light (4) with the on/off switch (8) on the base of the microscope.

**Experiments**

Use the following web link to find interesting experiments you can try out.

<http://www.bresser.de/downloads>

### TIPS on cleaning

Remove the device from its energy source before cleaning (remove plug from socket

/ remove batteries).

Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

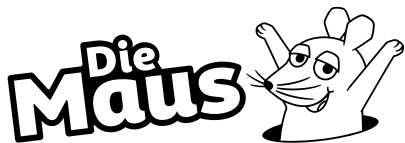
Clean the lens (objective and eyepiece) only with the cloth supplied or some other soft lint-free cloth (e.g. micro-fibre). Do not use excessive pressure - this may scratch the lens.

Dampen the cleaning cloth with a spectacle cleaning fluid and use it on very dirty lenses.

### Warranty & Service

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at [www.bresser.de/warranty\\_terms](http://www.bresser.de/warranty_terms).

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© I. Schmitt-Menzel
WDR mediagroup GmbH