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TABLE OF CONTENTS

IMPRINT	16
VALIDITY NOTE	16
ABOUT THESE INSTRUCTION MANUAL	16
1. GENERAL WARNINGS	16
2. SCOPE OF DELIVERY	17
3. CHOICE OF LOCATION & SETUP	18
4. PRODUCT DESIGN ADVANCE ICD 10X-160X ZOOM-STEREOMICROSCOPE	19
5. MICROSCOPE SETTINGS	21
6. CLEANING/MAINTENANCE/TRANSPORT	25
9. WARRANTY.....	26
10. CE DECLARATION OF CONFORMITY.....	26

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In the event of any warranty claims or service enquiries, please refer to the "Warranty" and "Service" sections of these operating instructions.

Subject to technical modifications without notice.

© 2024 Bresser GmbH. All rights reserved.

Reproduction of these operating instructions - in whole or in part - in any form (e.g. photocopying, printing, etc.) as well as use and distribution by means of electronic systems (e.g. image file, website, etc.) requires the prior written consent of the manufacturer and is otherwise prohibited.

The designations and brand names of the respective companies used in these instruction manual are protected by commercial, trademark and/or patent law in Germany, the European Union and/or other countries.

VALIDITY NOTE

These instruction manual is valid for products with the following article numbers:

5804000

MANUAL VERSION: 0424

MANUAL NAME:

Manual_5804000_Advance-ICD_de-en_BRESSER_v052024a

If you have any queries, please quote this information.

ABOUT THESE INSTRUCTION MANUAL

NOTE: These instruction manual are to be regarded as part of the device.

Read the safety instructions and the instruction manual carefully before using this device. Keep these instruction manual in a safe place for future reference. If the device is sold or passed on, the instruction manual must be handed over to the new owner/user of the product.

1. GENERAL WARNINGS

ATTENTION!

This device is not intended for use by persons (including children) with impaired physical, sensory or mental abilities, or for lack of experience and/or knowledge, unless they are supervised by or have received instructions from a person responsible for their safety on how to use this device.

DANGER OF AN ELECTRIC SHOCK!

This device contains electronic parts that are powered by a power source (mains connection). Never leave children unattended when handling the device! Only use the device as described in the instruction manual, otherwise there is a RISK of an ELECTRIC SHOCK!

- Disconnect the power supply when you are not using the device, during long breaks, and before all maintenance and cleaning work (disconnect the mains plug).
- Place the device so that it can be disconnected from the power supply at any time. The power outlet should always be near the device and should be easily accessible, as the plug of the power cord serves as a part to disconnect the device from the mains supply.
- To disconnect the device from the mains, always pull the mains plug and never the cable!
- Check the device, cables and connections for damage before use.
- Do not use the device if is damaged or contains damaged live parts! Damaged parts must be replaced immediately by an authorized service centre.
- Only operate the device in a completely dry environment and do not touch the device with wet or damp body parts.

DANGER OF SUFFOCATION!

Improper use can lead to a RISK OF SUFFOCATION, especially for children! It is therefore imperative that you observe the following safety information.

- Keep packaging materials (plastic bags, rubber bands, etc.) away from children!
- This product contains small parts that can be swallowed by children!
- If small parts are swallowed, seek medical advice immediately!

FIRE/EXPLOSION HAZARD!

Improper use of this product poses a RISK OF FIRE/EXPLOSION! Always observe the following safety instructions to avoid fires and/or explosions.

- Do not expose the device to high temperatures. Use only the supplied AC adapter. Do not short-circuit the device or dispose of it in fire! Excessive heat and improper handling can cause short circuits, fires and even explosions!

⚠ RISK OF INJURY!

This device contains components and/or accessories that can cause minor or serious body injury if used improperly. It is therefore essential that you observe the following safety information in order to prevent body injuries.

- This device often requires the use of sharp-edged and pointed tools. For this reason, keep this device together with all accessories and tools in a place inaccessible to children!
- Keep chemicals and liquids out of reach of children! Don't drink! Clean hands thoroughly under running water after use. In case of accidental contact with eyes or mouth, rinse with water. In case of complaints, consult a doctor immediately and show the substances.

! DANGER OF PROPERTY DAMAGE!

Improper handling can cause DAMAGE to the appliance and/or the accessories! Therefore, use the device only in accordance with the following safety information.

- Do not disassemble the device! In case of a defect, contact your dealer. The dealer will contact the service centre and send the device for repair if necessary.
- Do not expose the device to high temperatures and protect it from water and high humidity!
- Do not subject the device to shocks!
- Only use accessories and spare parts that comply with the technical specifications!

2. SCOPE OF DELIVERY

- A: Zoom stereo microscope with built-in LED incident and transmitted light illumination
 B: 2 pairs of wide-field eyepieces (WF 10x / WF 20x)
 C: 1 pc. Attachment lens (additional lens) (2x)
 D: 1 pair of rubber eyecups
 E: Plastic object plate, 94 mm (black/white)

- F: Object plate glass, 94 mm (matte)
 G: Power adapter (5V,1A) and 2 pcs. Plug adapter (1x EU, 1x UK) for the power adapter
 H: USB power cable (1 m)
 I: Dust cover
 J: Tool wrench for adjusting the friction

2.1. CHECK SCOPE OF DELIVERY / HANDLING TIPS

- Open the packaging with care. Avoid fingerprints and sweat on the optical surfaces.
- Prevent the microscope and accessories from falling and being damaged.
- Remove all components from the packaging and check for completeness according to the delivery note. Depending on the model or accessories, the microscope may be packed in more than one box.
- Handle the microscope and its accessories with care and protect them from strong impacts and vibrations at all times.
- Keep the original packaging for storage purposes or returning the microscope to the supplier in case repair or inspection is required.

ATTENTION! Never lift the microscope by the trinocular microscope head or the coarse/fine focusing knobs. This can damage the microscope.

ATTENTION! Always carry the microscope with both hands. To do this, hold the microscope with one hand on the underside of the microscope base while gripping the stand with the other hand.

3. CHOICE OF LOCATION & SETUP

Before setting up your microscope, choose a suitable location.

Place the microscope on a surface that meets the following criteria:

- Flat/level
- Vibration-free/stable
- Hard, non-combustible surface
- Chemically and mechanically resistant

Set up the microscope so that ...

- ... there is a distance of at least 10 cm around the device or to other appliances or the wall.
- ... you can disconnect it from the power supply at any time.
- ... The mains cable is visible and freely accessible at all times. The plug of the mains cable serves as a disconnecting device.
- ... the residual moisture has been completely removed before the first application.

ATTENTION! Do not expose the device to temperatures above 40° C! This can lead to irreparable damage to optical and/or mechanical parts that are not covered by the warranty.

ATTENTION! Always use the supplied power adapter (5V/1A) for the LED lighting. Using an unsuitable power supply unit can cause damage to the microscope!

4. PRODUCT DESIGN ADVANCE ICD 10X-160X ZOOM-STEREOMICROSCOPE

4.1. FRONT/SIDE VIEW

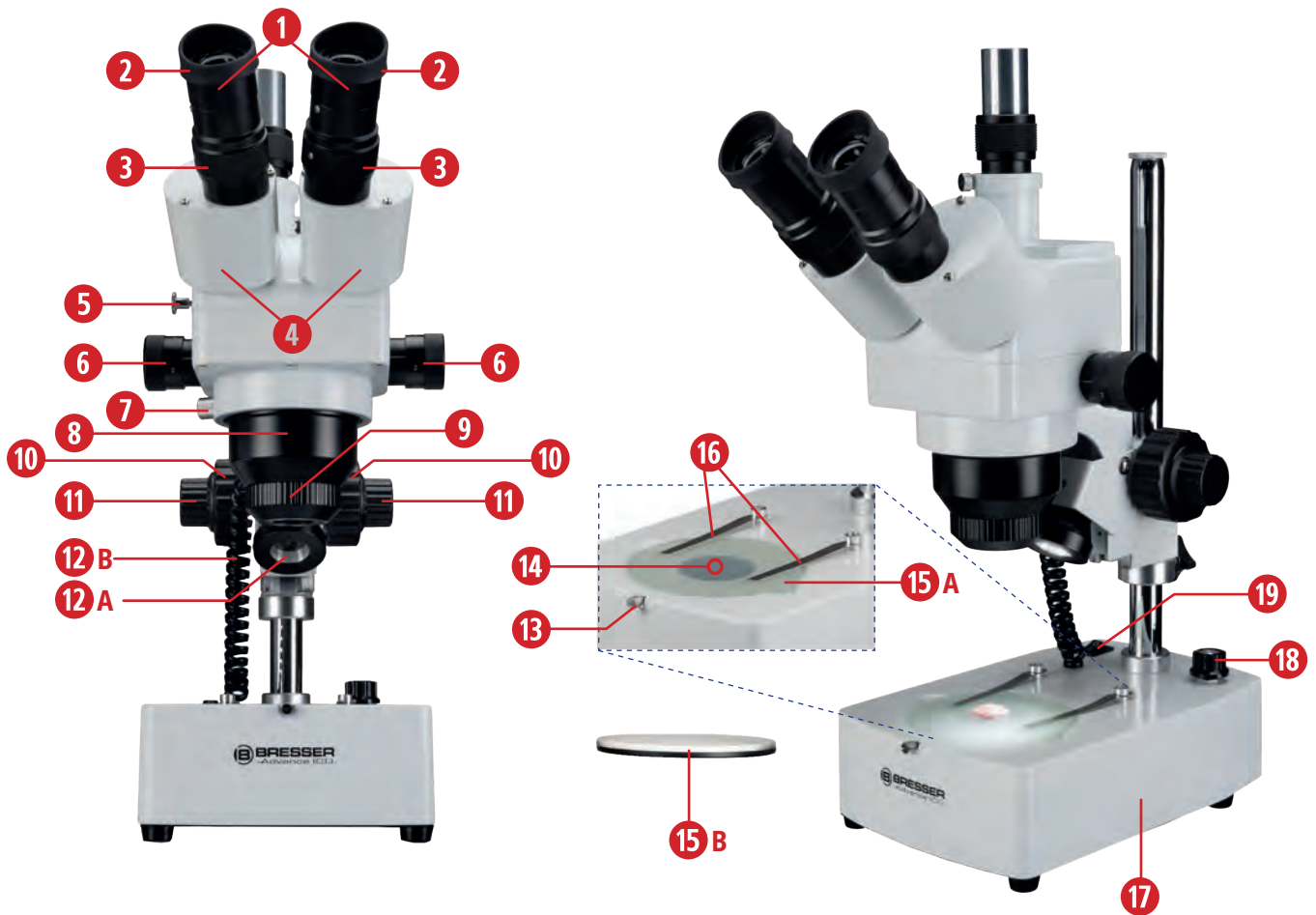
Fig. 1a: Components on the Advance ICD 10x-160x zoom stereo microscope front view.

EQUIPMENT

- | | |
|---|--|
| 1. Wide-field eyepieces (10x, 20x) | 12A. Reflected light unit |
| 2. Eyecups (rubber) | 12B. Connection cable for incident (reflected) light |
| 3. Diopter adjustment rings | 13. Locking screw-object plate |
| 4. Binocular attachment | 14. Transmitted light unit |
| 5. Lever bino/trino conversion | 15A. Glass object panel |
| 6. Zoom rotary knobs (1x to 4x) | 15B. Plastic object plate black/white |
| 7. Zoom lens retaining screw | 16. Clamp holder for preparation |
| 8. Stereo zoom lens (1x to 4x) | 17. Microscope base |
| 9. Additional lens (attachment lens) 2x | 18. Incident light/transmitted light switch |
| 10. Coarse focusing (coarse drive) | 19. On/Off switch |
| 11. Fine focusing (fine drive) | |

4.2. BACK VIEW

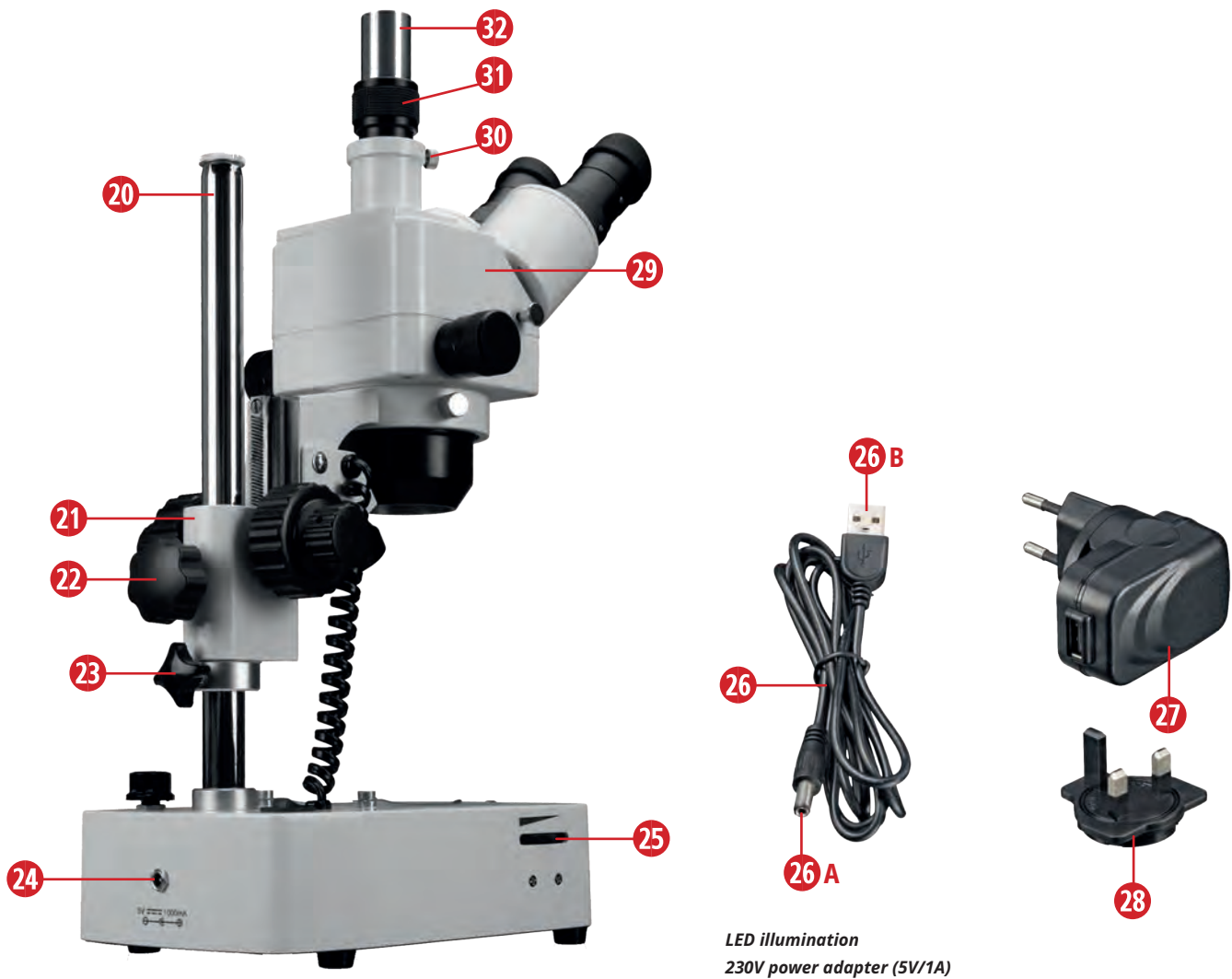
Fig. 1b: Components on the Advance ICD 10x-160x zoom stereo microscope Rear/side view.

EQUIPMENT

- 20. Tripod
- 21. Microscope arm
- 22. Microscope arm locking screw
- 23. Stopper with locking screw
- 24. mains connection socket
- 25. Dimmer for LED lighting
- 26. USB power cable with A) hollow plug and B) USB plug
- 27. Power adapter (5V/1A) with fitted EU plug adapter
- 28. UK adapter
- 29. Microscope head
- 30. Locking screw for adapter
- 31. Adapter with knurled ring
- 32. Trinotube for camera connection

5. MICROSCOPE SETTINGS

A POWER SUPPLY & ON/OFF SWITCH

I) Establish power supply

To establish the power supply, insert the hollow plug (1) of the supplied USB power cable (2) into the power connection socket (3) on the back of the microscope. Depending on the local area, you can fit the appropriate plug adapter (5) to the power adapter. Then plug the power adapter (4) into a suitable socket (7) (max. 230V).

Note: Two additional plug adapters (6) are included for the power adapter. Attach the plug adapter suitable for your socket to the power adapter (4).

ATTENTION! Make sure in advance that the plug adapter is firmly attached to the power adapter and cannot be detached from it! Never insert the plug adapter into the socket individually!

Fig. 2.: Establish power supply.

II) Switching the microscope on/off

To switch on the microscope, move the on/off switch (main switch) (4) to the 'I' position.

To switch the microscope off, move the ON/OFF switch to the '0' position.

B POSITION THE SAMPLE

I) Positioning of the sample

Place a sample (1) that you wish to observe on the glass plate (microscope stage) (2) and fix it in place with the clamps (3) if necessary.

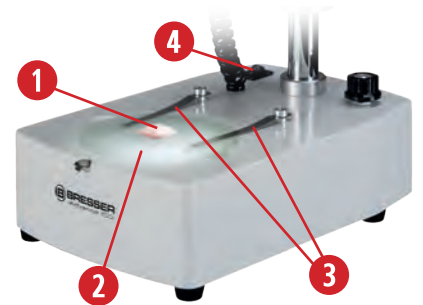
Fig. 3.: Position the specimen.

II) Use of the interchangeable object plates

The two object plates serve as a microscope stage. In addition to the frosted glass object plate (1), a further plastic object plate (2) with one black and one white side is also included in the scope of delivery. You can use this plastic plate to increase contrast depending on the object.

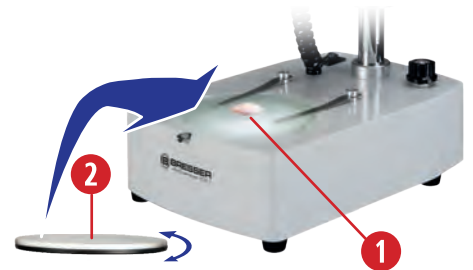
Fig. 4: Replacement of the object plate.

C SETTING THE LIGHTING

I) Switching the lighting on/off

You have the option of selecting three different illuminations for the object by setting the corresponding setting on the reflected light/transmitted light switch (1). In position "I", only the reflected light unit (2) is activated. The height of the incident light unit can be adjusted slightly to achieve the best possible illumination. In position "III", only the transmitted light unit (3) is switched on. In position "II", both lighting units are switched on. In the "OFF" position, the lighting is switched off.

II) Dimming the lighting

As the device is equipped with a continuously adjustable LED light (dimmer) (4), optimum illumination of the viewed object or sample is guaranteed.

ATTENTION! Do not leave the dimmer at maximum brightness for long periods of time. This shortens the lifespan of the lamp.

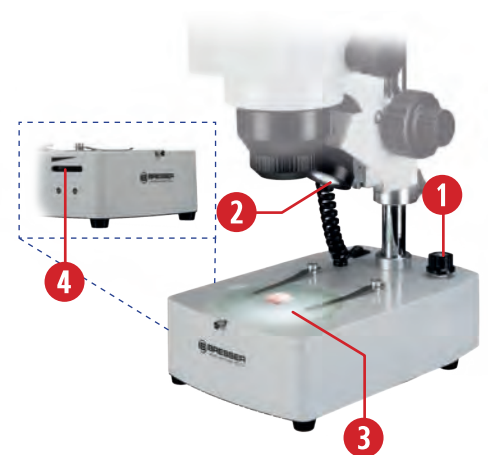
Fig. 5: Set the lighting type and dimmer.

Fig. 6: Adjust the interpupillary distance.

D ADJUSTING THE OBSERVATION TUBE

I) Adjusting the interpupillary distance (A) (distance between the eyepieces)

Set the correct interpupillary distance for binocular observation by shifting the eyepiece tubes (1) on the microscope head so that the right and left fields of view overlap completely or combine to form a circular image.

Remember: The correct pupil distance is crucial for comfortable, fatigue-free viewing.

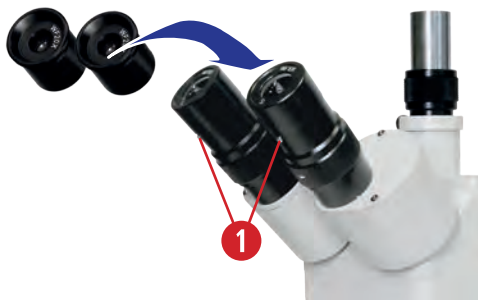
Fig. 7: Interchangeable eyepieces.

II) Replacing the eyepieces

You have two pairs of eyepieces available: WF 10x / WF 20x. You can easily replace these

To do this, loosen the eyepiece locking screws (1) with a small screwdriver. Make sure that you do not unscrew the screws too far. Once you have replaced the eyepieces, you can carefully tighten the screws again.

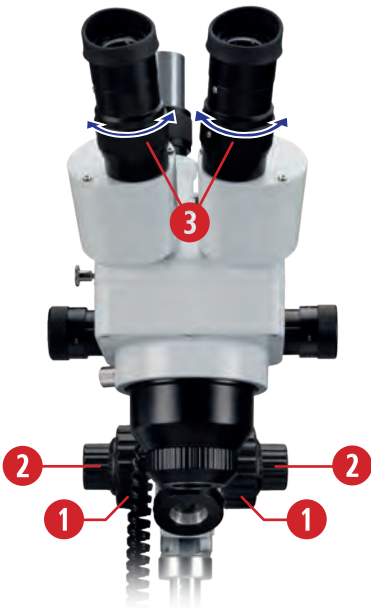
Fig. 8: Setting the diopter adjustment.

III) Adjusting the visual acuity correction (diopter compensation)

Focus a specimen as flat as possible using the coarse (1) and fine drive (2).

Now correct the sharpness of the left eye by closing the right eye and adjusting the diopter adjustment ring (3) of the left eyepiece accordingly until the left eye sees the sample sharply.

If necessary, proceed accordingly with the right eye.

Note: Each time you change the magnification using the zoom knobs, you adjust the focus using the focusing drives.

Note: Diopter compensation is carried out via the adjustment ring on both eyepieces (3) and is used to compensate for short-sightedness or far-sightedness (max. +/- 8 dpt. difference between the two eyes can be compensated). If the difference is higher or if astigmatism etc. is present, we recommend observation with your prescribed spectacles.

Note: diopter compensation is possible with both eyepieces, but is only performed for one eye if the difference is +/- 4 dpt. or less.

Fig. 9: Soft rubber eyecups.

IV) Use of rubber eyecups

The soft eyecups (1) included in the scope of delivery can be attached to the eyepieces. On the one hand, they protect spectacle wearers from damage in the event of accidental contact with the eyepiece and, on the other, they prevent distracting reflections on the lens surface caused by ambient light. They also keep the eyes at the correct distance from the eyepiece lens.

E FOCUSING THE SAMPLE

I) Set the image sharpness for a specimen (focus).

The microscope has a coaxial coarse and fine focusing mechanism. Turn the coarse drive (1) to raise or lower the microscope head (2). To make precise adjustments to the image sharpness, use the fine focusing knob (3).

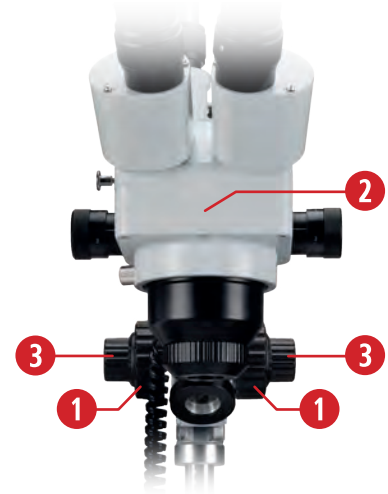
Fig. 10: Focusing the specimen.

II) Adjustment of the coarse drive (friction)

It can happen that the microscope head is no longer held by the focusing drive and moves down due to its own weight or heavy camera accessories. You can use the supplied tool wrench (1) to adjust the movement of the coarse drive (friction). To do this, select the coarse drive (2), which has an adjustment ring (3) with small round holes (4) on the inside. You can hook the tool wrench into these holes and turn the ring. Turning the ring towards the microscope makes the coarse drive run more stiffly and turning the ring away from the microscope makes the coarse drive run more smoothly.

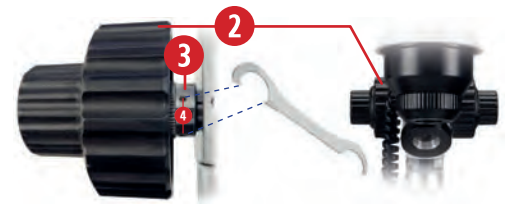
Fig. 11: Adjusting the movement of the coarse drive (friction).

III) Manual zoom (1x to 4x)

The stereo microscope has a stereo zoom lens (1). The magnification can be continuously adjusted between 1x and 4x by turning the zoom knobs (2).

Note: You can calculate the actual magnification by multiplying the value of the zoom by the value of the eyepiece (10x or 20x).

Note: Additional lenses change the working distance, whereby magnifications <math><1.0x</math> increase the working distance. This can be particularly advantageous when working under a microscope or handling tools. Magnification factors >math>>1.0</math> reduce the working distance.

Fig. 12: Stepless adjustment of the magnification.

IV) Use of the additional lens (attachment lens) 2x

You can additionally increase the magnification of your microscope by attaching the supplied 2x additional objective (1) to the stereo zoom objective (2). Carefully screw the additional lens into the mount provided on the lens.

Note: The combination of the stereo zoom objective, interchangeable wide-field eyepieces (10x, 20x) (3) and the 2x additional objective results in a very wide magnification range from 10x to 160x.

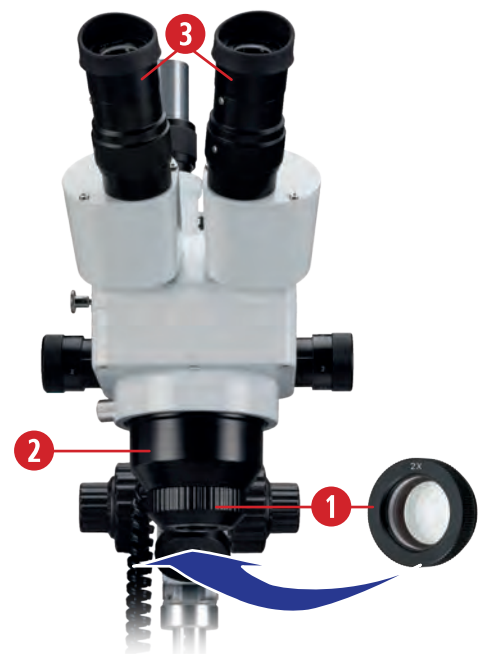
Fig. 13: Attaching the additional lens.

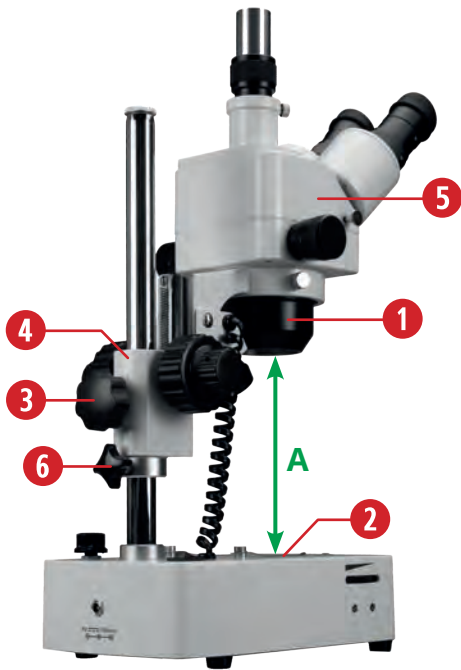
Fig. 14: Viewing of larger objects is possible via the height-adjustable microscope arm.

F VIEWING LARGER PREPARATIONS

The height-adjustable and 360° rotatable attachment also allows viewing of larger specimens, e.g. minerals or technical components, from any observation position.

The distance (A) between the zoom lens (1) and the object plate (microscope stage) (2) can be changed for this purpose. To do this, open the locking screw of the microscope arm (3) and adjust the distance according to the size of the object to be viewed. Then tighten the locking screw of the microscope arm again.

The microscope arm (4) holds the microscope head (5) and the zoom lens. The entire optical unit can be secured against sliding down using the locking screw of the stopper (6). Loosen the locking screw of the stopper and slide the stopper under the microscope arm. Then retighten the locking screw of the stopper.

NOTE: Maximum object height approx. 230 mm, working distance: 28-80 mm

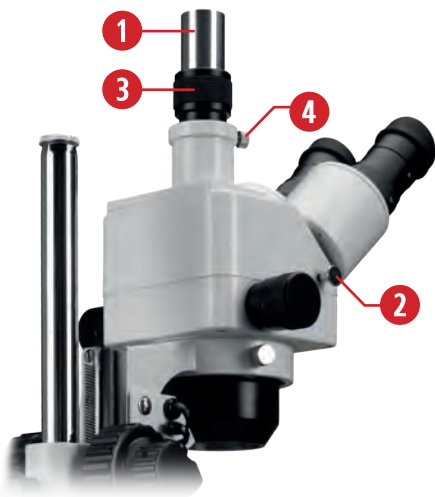
Fig. 15: Trinocular tube for photography

G TRINOCULAR TUBE FOR PHOTOGRAPHY

ATTENTION! Please find out in advance whether an additional adapter or a reduction lens is required for the use of the various cameras.

Use the height-adjustable trinotube (1) for photographic purposes. Various cameras can be connected. You can insert an eyepiece camera, such as the Bresser MikrOkular, directly into the trinotube without an additional adapter. The connection of a Bresser MikroCam may require additional adapters or reducing optics. To direct the beam path of the left eyepiece to the trinotube, pull out the bino/trinotube conversion lever (2) completely.

NOTE: The camera image may have a different degree of sharpness than the image produced at the eyepiece. Focus both images by adjusting the height of the trinotube using the knurled ring of the adapter (3).

Then fix the desired height of the trinotube using the adapter locking screw (4).

NOTE: The dust cap on the trinotube provides protection against dust and dirt. Remove the cap before photomicrography and reattach it after use.

6. CLEANING/MAINTENANCE/TRANSPORT

During long breaks, and before carrying out any maintenance and cleaning work, disconnect the power supply by unplugging the mains plug.

Remove all interchangeable optical elements (e.g. additional objectives, eyepieces, etc.) before cleaning.

As the microscope is a precision instrument, please do not dismantle any parts, otherwise proper use can no longer be guaranteed.

CLEANING

First, blow loose dust off the lens surfaces.

Use a soft disposable paper towel soaked in 30% to 70% pure alcohol (available from pharmacies or drugstores). Do not apply the cleaning solution directly to the optical parts. Gently wipe the lens surface by cleaning the surface with circular movements from the center to the edge of the lens. Apply light pressure to the optics.

ATTENTION! Do not use organic solvents (e.g. alcohol, ether, acetone, xylene or other thinners) to clean painted parts or plastic parts!

ATTENTION! Alcohol is highly flammable.

ATTENTION! Do not apply cleaning fluid directly to the appliance to avoid damaging the electronics.

Dust cover:

When not in use, cover the microscope with the dust cover and place it in a dry and mold-free place. We recommend storing all objectives and eyepieces in a closed container with desiccant.

Transport:

Remove all loose components and the specimen to be observed from the cross table. Please pack the microscope carefully (if possible in the original packaging).

7. TECHNICAL DATA

- Trinocular attachment for visual observation and simultaneous camera use
- Various cameras can be connected
- 1x to 4x stereo zoom lens, magnification: 10x-160x
- 2x additional lens (doubles the magnification with eyepieces and zoom lens)
- Bright LED lighting
- Two different object plates for optimum contrast to the object
- Observation of three-dimensional objects up to a maximum object height of approx. 230 mm possible
- Dimensions (LxWxH): 330 x 148 x 620 mm / weight: 4,7kg
- Power adapter 5V/1A, a matching USB power cable with a length of 1 m and two plug adapters (EU and UK) are included in the scope of delivery
- Solid design with heavy base
- 2 pairs of wide-field eyepieces (10x/20x) included

Eyepiece:	Magnification:	Field of view:
10x wide field eyepiece	10 - 40x	approx. 22 - 5.5 mm
10x wide field eyepiece with additional 2x lens	20 - 80x	approx. 11 - 2.5 mm
20x wide field eyepiece	20 - 80x	approx. 11 - 2.5 mm
20x wide field eyepiece with additional 2x lens	40 - 160x	approx. 5.5 - 1.1 mm

8. DISPOSAL

 Do not dispose of electrical appliances with household waste! In accordance with Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its transposition into German law, used electrical appliances must be collected separately and recycled in an environmentally friendly manner.

 Please observe the current legal regulations when disposing of the microscope and all accessories! Information on proper disposal can be obtained from municipal waste disposal service providers or the Environmental Agency.

Dispose of the packaging materials according to type. Information on proper disposal can be obtained from municipal waste disposal service providers or the Environmental Agency.

9. WARRANTY

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as indicated on the gift box, registration on our website is required.

You can view the full warranty conditions and information on warranty extensions and services at www.bresser.de/warranty_terms.

www.bresser.de/warranty_terms

10. CE DECLARATION OF CONFORMITY

CE Bresser GmbH has prepared a "CE Declaration of Conformity" in accordance with the applicable guidelines and corresponding standards. The full text of the CE Declaration of Conformity is available at the following website: www.bresser.de/download/5804000/CE/5804000_CE.pdf

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

E-Mail: service@bresseruk.com

Telephone*: +44 1342 837 098

BRESSER UK Ltd.

Suite 3G, Eden House

Enterprise Way

Edenbridge, Kent TN8 6HF

Great Britain

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Bresser UK Ltd.