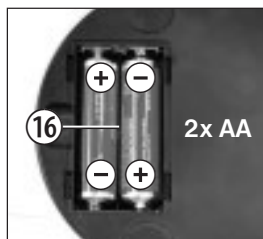
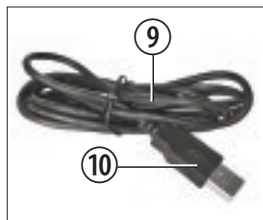
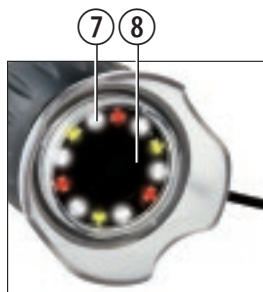
UNIVERSAL MIKROSKOP UNIVERSAL MICROSCOPE 20X/200X

- DE** Bedienungsanleitung
- EN** Operating instructions
- FR** Mode d'emploi
- NL** Handleiding
- IT** Istruzioni per l'uso
- ES** Instrucciones de uso
- RU** Руководство по эксплуатации

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Kaufdatum:

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General Warning

- This device contains electronic components that operate via a power source (power supply and/or batteries). Only use the device as described in the manual; otherwise, you run the risk of an electric shock.
- Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short circuit, a fire or an explosion.
- Never bend, pinch, pull or run over the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Children should only use the device under adult supervision. Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as they pose a choking hazard.
- Before cleaning the device, disconnect it from the power supply by removing the batteries.
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.

Declaration of Conformity

 Bresser GmbH has issued a 'Declaration of Conformity' in accordance with applicable guidelines and corresponding standards. This can be viewed any time upon request.

Parts

1. Camera shutter
2. Reduce brightness (-)
3. Increase brightness (+)
4. Set light mode/turn light off
5. Focus ring
6. Attachment piece
7. Illuminator (12 LEDs)
8. Lens
9. USB cable
10. USB connector
11. Stand
12. Focusing knob
13. Illuminated table
14. Preparation holders
15. Box with object holder,
glass covers and adhesive labels
16. Battery compartment (stand)

General

This is a digital reflected-light microscope. You hold it in your hand and can place the bottom section (attachment piece) on all kinds of things in order to look at them. For example, you can observe leaves, microorganisms, your skin or hair and much more. The microscope works best when the thing that you're looking at (also called the 'object') is flat. You can view the enlarged image on your computer screen, as well as take still pictures and video and save them on your computer.

Driver Installation

Insert the product CD into your PC's CD/DVD drive. The driver installation starts automatically. Plug the microscope's USB connector (10) into your computer's USB slot. The lighting (7) turns on and your PC detects the hardware, which is then installed. Soon, the 'AMCAP' icon appears on the desktop. Now you can use the microscope.

Live Observation

Press the camera shutter (1). A live image is displayed on the monitor. It will probably be blurry. Hold the microscope by the casing and place the attachment piece (6) on an object, like a piece of paper with writing on it. Turn the focus ring (5) to make the live picture sharper. This is called focusing.

Light Mode Settings

Press the 'MODE' key to select from four different color combinations:

- white light
- white combined with red light
- white combined with yellow light
- white combined red and yellow light

Thus you will receive the optimal illuminated light depending on the object. The light can be shut off altogether by pressing the 'MODE' key as well. Regulate the brightness using the '+' and '-' keys on the device by pressing and holding the keys.

Taking Pictures

Using the camera shutter (1), you can take a picture and save it as a BMP file.

1. Press the camera shutter (1).
2. The 'SnapShotView' window appears on the screen with a picture.
3. To save the image, click 'File' and 'Save'.

Making Movies

The 'AMCAP' program allows you to make movies with the microscope and save them as AVI files.

1. Click on 'File' and 'Set Capture File...'; specify the name of the AVI file with the file extension '.avi'. For example: 'experiment1.avi'.
2. The 'Set File Size' window appears on the screen. Here, specify the maximum file size.
3. You can prepare to film with 'Capture' and 'Start Capture' in the menu.
4. Start filming with 'OK' in the 'Ready to Capture' window.
5. Under the 'Capture' menu, you can end your filming with 'Stop Capture'.
6. If you want to record a new film, follow Step 1 and specify a new AVI file with a new name. Otherwise, the file will overwrite your film.
7. You can watch your film using a playback program for multimedia files.

Magnifications

In the lower magnification, a picture includes about 10.5 mm x 14 mm of the object. The higher magnification includes about 1 mm x 1.4 mm. The higher mag-

nification is about five times stronger than the lower one. When, for example, you print a picture that is 28 cm wide on a piece of paper, the magnification is about 20X (low) or 200X (high).

Deactivation and Storage

Close the 'AMCAP' window on your PC screen. Now, you can remove the USB connector (10) from your computer's USB port. You can store your microscope in the product packaging. This will protect it from dust.

Technical Information

- Digital handheld microscope with computer connection (USB)
- Magnification: 20X and 200X
- Bright illumination via 12 LEDs
- Power supply via USB
- Size: 54 x 54 x 104 mm
- Weight: 144 g

System Requirements

Windows XP with Service Pack 3 (on CD-ROM), Windows Vista, Windows 7 with DirectX 9.x (on CD-ROM), a minimum of 1 GB RAM, free USB 2.0 port

Photomizer SE Software

Photomizer SE Software can be downloaded free of charge from <http://www.bresser.de/downloads/support/software/photomizer.zip>

Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

Empty, old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 6 January 2006 from your local waste-disposal service or environmental authority.

 In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law – at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

Batteries that contain toxins are marked with a sign and a chemical symbol.

Cd¹

Hg²

Pb³

¹ battery contains cadmium

² battery contains mercury

³ battery contains lead

Warranty and warranty term extension

The warranty term is two years from the date of purchase. Please keep your proof of purchase. Register at www.bresser.de/warranty and fill out a brief questionnaire to get your warranty term extended to **five years**. Registration must be completed within three months of purchase (date of receipt) to validate the warranty. If you register thereafter, the warranty term will not be extended.

If you have problems with your device, please contact our customer service first. Do not send any products without consulting us first by telephone. Many problems with your device can be solved over the phone. If the problem cannot be resolved by phone, we will take care of transporting your device to be repaired. If the problem occurred after the warranty ended or it is not covered by our warranty terms, you will receive an estimate of repair costs, free of charge.

Service Hotline: +49 (0) 2872 - 80 74-210

Important for any returns:

Please make sure to return the device carefully packed in the original packaging in order to prevent damage during transport. Also, please enclose your receipt for the device (or a copy) and a description of the defect. This warranty does not imply any restriction of your statutory rights.

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are trademarks of the National Geographic Society,
used under license. All rights reserved.