- DE** Bedienungsanleitung
- EN** Instruction Manual
- FR** Mode d'emploi
- ES** Instrucciones de uso

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General Information

About this Instruction Manual

Please read the safety instructions in this manual carefully. To avoid damage to the unit and the risk of injury, only use this product as described in the manual. Keep the instruction manual handy so that you can easily look up information on all the functions.

**DANGER!**

You will find this symbol in front of every section of text which deals with the risk of severe injury or even death in the event of improper use.

**CAREFUL!**

You will find this symbol before every section of text that deals with the risk of minor to severe injuries resulting from improper use.

**NOTE!**

You will find this symbol in front of every section of text which deals with the risk of damage to property or the environment.

**CAREFUL!**

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after use. In the event of contact with the eyes or mouth, rinse thoroughly with water. In the event of pain, contact a doctor immediately and take the substances with you.

**NOTE!**

Do not disassemble the device. In the event of a defect, please contact the Service Centre in your country (see the warranty card).

Do not expose the device to temperatures above 40°C.

Intended Use

This product is intended only for private use. It was developed for the magnified display of things in nature.

**General Warning****DANGER!**

Tools with sharp edges and points are often used when working with this device. For this reason, store this device and all accessories and tools in a location that is out of the reach of children. There is a RISK OF INJURY!

**DANGER!**

This device contains electronic components which operate via a power source (power supply and/or batteries). Do not leave children unattended when using the device. Only use the device as described in the manual, otherwise you run the RISK of an ELECTRIC SHOCK.

**DANGER!**

Do not expose the device to high temperatures. Use only the supplied power adapter. Do not short circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short circuit, fires or even explosions.

**DANGER!**

Never bend, pinch, pull or run over the power and connecting cables or extensions/adapters. Protect the cables from sharp edges and heat. Before operation, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables, etc. Damaged parts must be exchanged by an authorized service centre immediately.

**DANGER!**

Children should only use the device under adult supervision. Keep packaging material (plastic bags, rubber bands, etc.) out of the reach of children! There is a RISK OF CHOKING!

1. Components

- 1 Eyepiece
- 2 Eyepiece holder
- 3 Microscope head
- 4 Objective revolver
- 5 Objective
- 6 Slide clamp
- 7 Condenser
- 8 Iris diaphragm (with adjusting lever)
- 9 Filter holder
- 10 Condenser height adjustment knob
- 11 Illumination
- 12 Coarse-adjusting wheel
- 13 Fine-adjusting wheel
- 14 Lengthwise knob of mechanical desk
- 15 Cross knob of mechanical desk
- 16 Dimmer
- 17 On/Off switch
- 18 Status-LED (red, green)
- 19 AC power socket

2. Package contents

(several accessory parts are premounted)
Depending on the version a 60x (5102060) or 100x (5102000) objective is included.

Microscope head
10x WF eyepiece
Objectives: 4x, 10x, 40x, 60x or 100x (oil)
Condenser
Filter plate (blue)
AC adapter
Dust proof cover

3. General/Location

Before you set up the microscope, you must choose a suitable location.

First you must make sure that your microscope is on a stable and solid surface.

4. Electric illumination

It works, by plugging the connection cable into the microscope socket (19) and plugging the AC adapter into the mains socket. The switch at the rear of the microscope must be switched on. HINT for the rechargeable battery:

Your microscope is fitted with one built-in rechargeable battery. A fully charged battery allows ca. 48 hours of observation. Before using the microscope for the first time the rechargeable battery should be charged for 16 hours. Please charge it by connecting the microscope socket (19) with the included AC adapter and plugging the AC adapter into mains socket.

5. Observation

You begin each observation with the lowest magnification (eyepiece 10x and objective 4x). So, the object to be viewed can easily be centered and properly focussed. The higher the magnification the more light is required for good image quality. For applications with the 100x lens, it is recommended that examinations are done only with oil immersion.

Begin with a simple observation. Turn the objective revolver (4) onto the smallest magnification and insert the 10x eyepiece. Now place a preparation directly under the objective on the microscope table. Move the preparation on the table with the two adjusting screws (14 and 15).

Now look through the eyepiece (1) and turn the coarse adjustment (12) carefully until the image becomes visible. The optimal sharpness is achieved by the fine adjustment (13).

Make sure that you never overwind the fine adjustment.

6. Condenser

Use the condenser height adjustment knob (10) to properly concentrate the light for your object.

7. Iris diaphragm

An iris diaphragm (8) is available if you want better image quality. This makes it possible to change perspective depth. The smaller the aperture the greater the perspective depth. However, image brightness is also reduced. You can compensate for this by adjusting the brightness using the dimmer (16).

8. Filter holder

Push the filter holder (9) forwards and then insert the blue filter if you wish to use it. The filter holder can then be pushed back again. This filter is used to show natural colours.

9. Oil immersion (only for 100x objective)

The immersion oil is only necessary for the 100x objective. Apply one drop of immersion oil onto the upper condenser lens. With the slide clamp (6), fix your prepared slide on the mechanical desk. Using the condenser height adjustment knob (10), move up the condenser (7) carefully, until the slide touches the immersion oil. Focus the preparation with the 100x objective and the fine focuser as good as possible. Now turn the objective revolver a little bit to the side and apply one drop of immersion oil onto the preparation's covering glass. Turn back the objective revolver. The objective must dip into the oil! Now the preparation can be focused. Don't forget to clean objective and preparation with alcohol after observing.

NOTES on cleaning

1. Avoid disassembling the microscope!
Because it is a precise optical instrument, do not disassemble the microscope casually. That may cause serious damage to its performance. Do not disassemble the objectives.
2. Avoid contamination!
You should make sure that dust or moisture do not get in contact with your microscope. Avoid touching optical surfaces with your fingers! If dust or dirt nevertheless have gotten onto your microscope, or the accessories, remove it as described below.
3. Cleaning non-optical parts:
Disconnect the plug from mains socket before cleaning non-optical parts of the microscope (e. g. the body or the stage). First, remove dust or dirt from your microscope, or accessories, using a soft brush. Then clean the dirty area with a soft, lint-free cloth; the cloth may be moistened with a mild detergent solution. Be sure that the microscope is dry before using. For cleaning the painted parts or plastic components, do NOT use organic solvents such as alcohol, ether, acetone, xylene or other thinner etc.!
4. Cleaning optical parts:
In order to enhance the optical quality, the lenses of eyepieces or objectives are coated. They should not be wiped because dry dirt or dust may scratch the coating. It is best to remove the parts to be cleaned from the body prior to cleaning. Always blow loose dust away first. Use lens tissue of good quality or soft fabric moistened with a small amount of pure alcohol (available in the pharmacy); wipe the lens surface clean.
5. Cleaning the oil immersion lens:
After working with oil immersion, the immersion oil should be removed from the lens of the 100x objective at the end of each workday. In order to do this, use a lens tissue moistened with a small amount of pure alcohol (available in the pharmacy). Don't forget to clean the preparation, too.
6. Storing the microscope:
After usage, cover the microscope with the dust cap (made of plastic) and keep it at a dry place, away from mold. If it is not used for a prolonged time, store the microscope and accessories in their respective containers.

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

Technical data

Monocular head
Mechanical desk (with vernier alignment)
Eyepiece: DIN WF 10x
Objectives: DIN 4x / 10x / 40x / 60x or 100x (oil)
Magnification: 40x - 600x or 1000x
Illumination: rechargeable battery LED

Technical alterations and errors excepted.

Aperture:

The numerical aperture (the unit of measurement for the resolution factor of the objectives) is engraved beside the magnification-sign on the objectives.

The magnification for microscopes

is calculated by formula:

Eyepiece magnification x objective magnification factor
= total magnification

Example: 10x eyepiece x 100x lens
= 1000x total magnification

Disposal

Dispose of the packaging materials properly, according to their type (paper, cardboard, etc). Contact your local waste disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage! As per the Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner. Empty old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 01.06.2006 from your local waste disposal service or environmental authority.

Please take the current legal regulations into account when disposing of your device. You can get more information on the proper disposal from your local waste disposal service or environmental authority.

Declaration of Conformity

Product Type: Transmission Microscope

Product Name: Erudit DLX

Article No.: 5102000 / 5102060

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address:

www.bresser.de/download/5102000/CE/5102000_CE.pdf

www.bresser.de/download/5102060/CE/5102060_CE.pdf

Bresser GmbH

DE-46414 Rhede/Westf. · Germany

Warranty & Service

The regular guarantee period is 5 years and begins on the day of purchase. You can consult the full guarantee terms and details of our services at www.bresser.de/warranty_terms.

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